CMP-USBMICRO10

MANUAL (p. 2)
USB MICROSCOPE

ANLEITUNG (s. 4)
USB-MIKROSKOP

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ENGLISH

How To Use The Holder

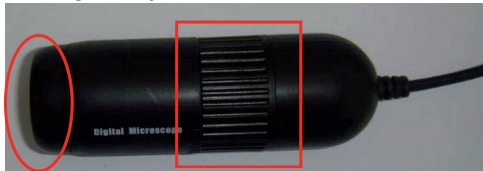
To change the angle of the holder, loosen the bolt carefully and move the holder gently. The holder may be broke when you try to move it while the bolt is tightened. The complete stand and holder is shown below.

Focus Adjustment

To watch a clear video image, be sure to rotate the focus ring to the right position to get a clear view. In the figure, a red box shows the focus ring of the microscope camera. The lens side of the microscope camera is indicated with a red oval. This end should be close to the object for getting a large magnification view.

Observing Correctly With The Stand


The stand is used to help you hold the microscope camera steady when observing an object. The common way to insert the microscope camera in the holder is shown in the picture below.

Software Installation

- Insert the CD-ROM or click on the icon of the CD device, then select "AutoPlay" option.
- Click "Next"
- Choose "I accept the terms of the license agreement" and Click "Next"
- Enter a user name and/or company name and product serial number, then click "Next".
The product serial is on the CD jacket.
- Click "Next". Note: Do not change the file installation folder.
- The program will install the files automatically.

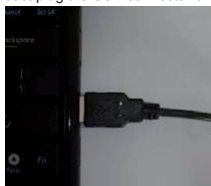- After the software installation is complete, an icon is added to your desktop. Double-click the icon to run the software for the microscope camera.
- Because this product uses an universal video driver, it is not necessary to install the camera driver. Just plug the USB connector of the microscope camera into a computer, as shown below.

A message will come up at the lower right corner of the computer screen, as below:
They show a microscope has been connected correctly and can be used.

For detailed information about the product and software, please check the full manual on the CD!

Safety precautions:

CAUTION
RISK OF ELECTRIC SHOCK
DO NOT OPEN

To reduce risk of electric shock, this product should **ONLY** be opened by an authorized technician when service is required. Disconnect the product from mains and other equipment if a problem should occur. Do not expose the product to water or moisture.

Maintenance:

Clean only with a dry cloth. Do not use cleaning solvents or abrasives.

Warranty:

No guarantee or liability can be accepted for any changes and modifications of the product or damage caused due to incorrect use of this product.

General:

Designs and specifications are subject to change without notice.

All logos brands and product names are trademarks or registered trademarks of their respective holders and are hereby recognized as such.

Keep this manual for future reference.

Attention:

This product is marked with this symbol. It means that used electrical and electronic products should not be mixed with general household waste. There is a separate collections system for these products.

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Copyright ©
NEDIS B.V.